


## TriPlus Crossrail X Installation Alignment

Only for Single TRACE GC Ultra and Single FOCUS GC

Loosen the fixing screws present on each support leg. Standing in front of the crossrail X let it slide gently until the marker present towards the crossrail left end is perfectly aligned with the reference marker placed on the leg of the left support.


The alignment of the relevant marker is recommended to allow the operator to take advantage of the predefined centering of sampling injection device on the injectors. The operator shall only perform the fine adjustment.

In all the other TriPlus-GC configurations the sampler injection device predefined centering is not available.

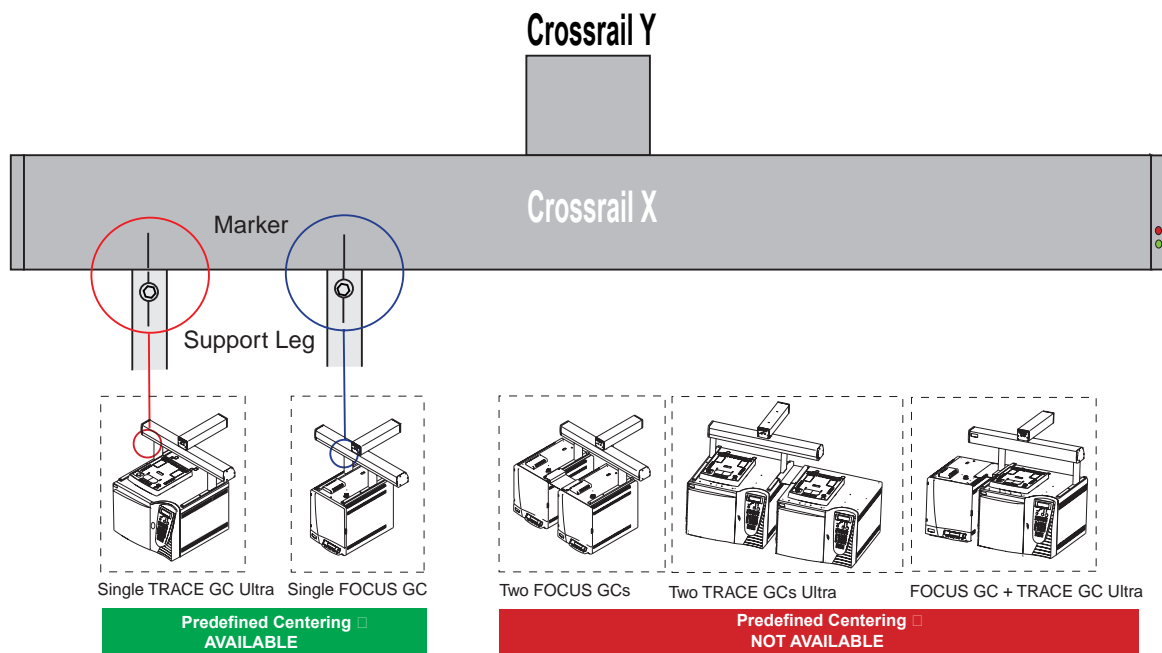

# TriPlus Components Installation Limits

The tables report the left and right installation limits of the components on the TriPlus Sampler. Each limit is measured from the left/right outer end of the crossrail X (standard / extended) and the left/right outer end of the component of interest.

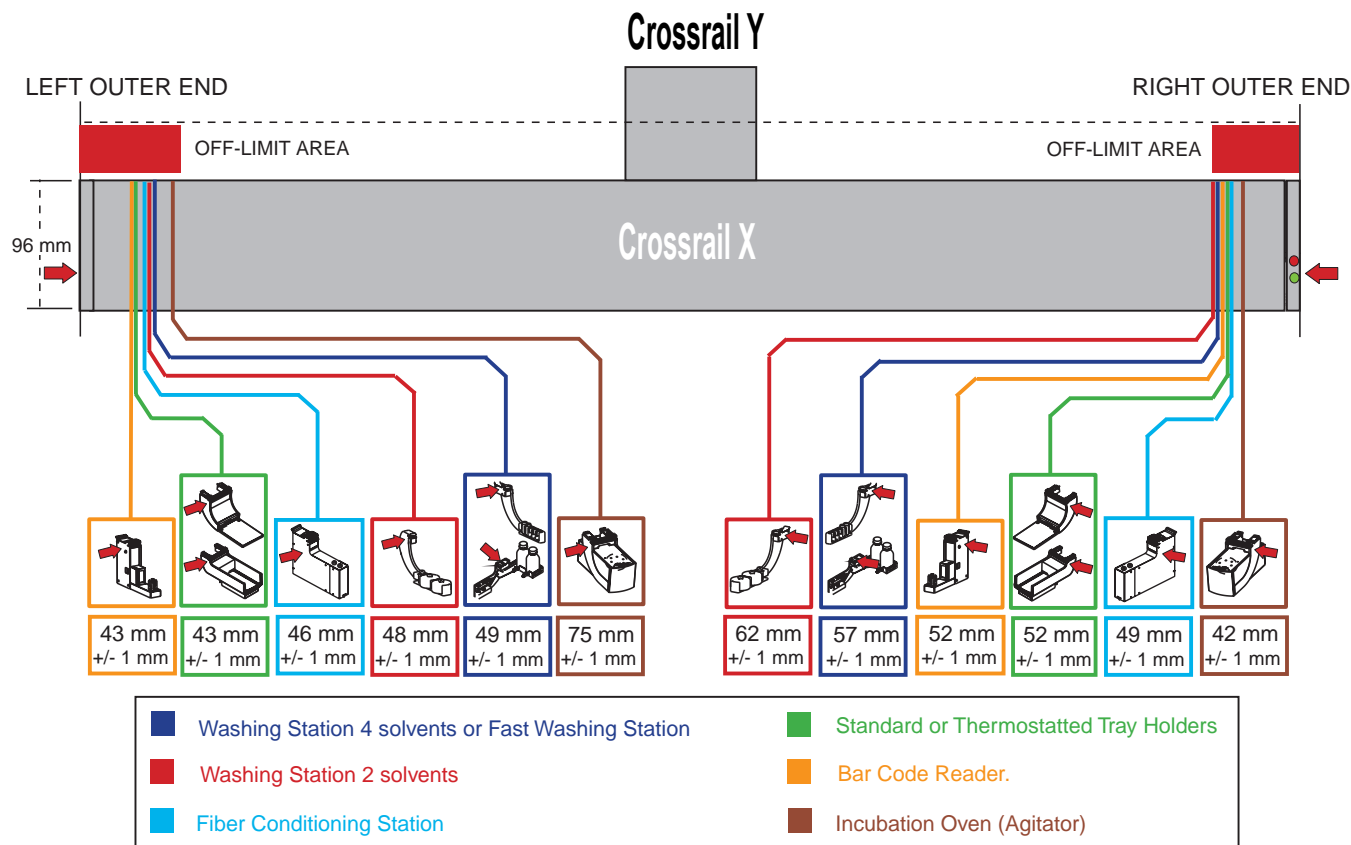

**CAUTION** The washing station module, when used in combination with **HS/SPME** version, requires the optional Vial bracket holder. The washing station equipped with Vial bracket holder is required to be mounted on the **LEFT SIDE** of the longitudinal crossrail of the sampler. The Vial bracket holder is not required when operating with **AS** version, and the washing station can be mounted along the longitudinal crossrail also on **RIGHT SIDE**.